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8-2 Error Code Table

Error code	Cause	Countermeasures
Er1	The sterilizing water is insufficient.	Add tap water.
	The water level sensor is contaminated.	Clean the water level sensor.
	Pure water is used for sterilizing water.	Use tap water for sterilizing water.
	The temperature of the sterilizing water is extremely low.	Press the CHECK key for approximately 30 seconds to increase the water temperature.
Er2	The exhaust valve is open.	Turn the exhaust valve knob clockwise to close the valve.
Er3	The temperature of the outer wall of the chamber is abnormally high during sterilizing.	Unplug the power cord and contact your dealer or the nearest TOMY office.
Er4	The volume of chamber water becomes insufficient during sterilizing.	Add sterilizing water after the chamber pressure falls 0kPa and the chamber temperature is 97°C or less.
Er5	The chamber temperature rises at least 5°C higher than the setting value during sterilizing.	Unplug the power cord and contact your dealer or the nearest TOMY office.
Er6	The pressure rises abnormally as the form of materials to be sterilized prevents air from escaping.	Turn off the power switch temporarily and wait for the chamber pressure to return to 0kPa and open the chamber lid to release air from the sterilizing material or to reduce the volume of the material. (See 5-6. Replacing the Material to Be Sterilized.)
Er7	The temperature sensor has fractures.	Unplug the power cord and contact your dealer or the nearest TOMY office.
Er8	The volume of the material to be sterilized is too much.	Open the chamber lid and reduce the volume of the material after the chamber pressure returns to 0kPa.
	The heater has fractures.	Unplug the power cord and contact your dealer or the nearest TOMY office.
Er9	The chamber lid is not securely seated in the chamber arm guide.	Reopen the lid and then close it firmly in place.
Er0	There are several potential causes stated above for the error.	Identify causes applying to the error referring to the above stated error codes from Er 1 to 9, and take appropriate actions corresponding to the relevant error codes.